

Thermosel™

ACCESSORY

Operator Instructions

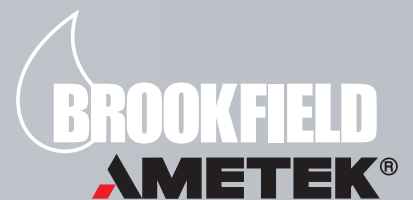


TABLE OF CONTENTS

I. Introduction	5
I.1 Utilities	5
I.2 Specifications	5
I.3 Safety Symbols and Precautions.....	6
I.4 Operation	6
I.5 General.....	10
I.6 Viscosity Measurements	10
I.7 Clean-Up	13
Appendix A - Viscosity Ranges / Spindle Factors	14
Appendix B - High Temperature Calibration Fluids	18
Appendix C - Disposable Sample Chamber And Spindles	19
Appendix D - Solid Shaft Spindle Option.....	21
Appendix E - Online Help and Additional Resources	22
Appendix F - Warranty Repair and Service	23

Thermosel™ Accessory Operator's Instructions

NOTE: These instructions apply to both the Standard and Magnetic versions.

I. INTRODUCTION

The Brookfield Thermosel System consists of a Brookfield Viscometer or Rheometer with related accessories to accurately measure the viscosity of liquids at elevated temperatures. Viscosity measurements are obtained using coaxial cylinder geometry. Temperature is controlled by a solid state, triac output, Proportioning Controller which maintains the spindle, chamber, and sample material at the desired temperature.

The system is designed for measuring liquid viscosities over a temperature range from 40°C (104°F) to 300°C (572°F). The designed viscosity range is from 1.2 cP to 80,000,000 cP depending upon the viscometer and SC4 spindle used.

Please note that all references made herein to digital viscometers would also apply to DV3T Rheometers and DVNext.

I.1 UTILITIES

Input Voltage:	115 VAC or 230 VAC
Input Frequency:	50/60 Hz
Current Rating:	15 Amps



Main supply voltage fluctuations are not to exceed $\pm 10\%$ of the nominal supply voltage.

I.2 SPECIFICATIONS

Operating Environment: 5°C to 40°C temperature range (32°F to 104°F)

Relative Humidity: 20% - 80% R.H.: non-condensing atmosphere

Conforms to CE Standards:

BSEN 61326-1:2013:	Electrical equipment for measurement, control and laboratory use - EMC requirements.
BSEN 61010-1:2010:	Safety requirements for electrical equipment, for measurement, control and laboratory use.
BSEN IEC 63000:2018:	Technical documentation for the assessment of electrical and electronic products with respect to the restriction of hazardous substances (ROHS 3).

I.3 SAFETY SYMBOLS AND PRECAUTIONS

Safety Symbols:

The following explains safety symbols which may be found in this operating manual.



(Hot Surface Symbol)

Indicates surface may be hot.



(Caution Symbol)

Refer to the manual for specific warning or caution information to avoid personal injury or damage to the instrument.

Precautions:



(Caution Symbol)

If this instrument is used in a manner not specified by the manufacturer, the protection provided by the instrument may be impaired.



(Caution Symbol)

This instrument is not intended for use in a potentially hazardous environment.



(Caution Symbol)

In case of emergency, turn off the instrument and then disconnect the electrical cord from the wall outlet.



(Caution Symbol)

The user should ensure that the substances placed under test do not release poisonous, toxic or flammable gases at the temperatures which they are subjected to during the testing.

I.4 OPERATION

VISCOMETER AND STAND

The instrumentation should be placed on a firm level surface near a 15 amp, 115 or 230 volt, 50 or 60 Hz grounded AC electrical service, as required. A cooling water supply and drain arrangement is also recommended for cooling the Thermo Container with optional cooling plug assembly.

Set up the viscometer stand by connecting the upright rod with clamp attached into the base. Screw the leveling screws into position on the base. (Refer to the assembly instruction supplied with your lab stand.) Attach the viscometer to the stand by inserting it into the clamp; level the viscometer and position it centered between the stand legs. Raise the viscometer on the lab stand upright rod so that there will be sufficient room to attach the alignment bracket. Check that the power switch is OFF and plug the viscometer power cord into the appropriate electrical service.

ALIGNMENT BRACKET

Attach the alignment bracket to the rear of the viscometer pivot cup, securing it tight with the knurled screw. (Refer to Figure I-2).

THERMO-CONTAINER, SAMPLE CHAMBER, SAFETY GUARD

Using the three (3) leveling screws in the Thermo Container base, level the base.

Check that the perforated safety guard is flush onto the base of the Thermo Container.

Using the extracting tool, insert the chamber into the thermo-container. Rotate the chamber until it drops and locks in place prohibiting further rotation.

CONTROLLER AND PROBE

Set the controller on a level surface adjacent to the Thermo Container. Insert the male plug (three-prong) from the Thermo Container braided cord into the socket on the back of the controller (turn and lock connection). Insert the four-inch stainless steel probe (Resistance Thermometer) into the hole in the Thermo Container located directly above the braided cord. Plug the other end of the probe into the connector located on the back of the controller. To remove the probe, depress the spring clip and carefully slide probe out. This removal procedure should not be followed if the Thermo Container is at a temperature above 100°F (37°C). Cool the Thermo Container first.



If the probe is not inserted in the Thermo Container and the controller is turned on, the Thermo Container will cease to heat and the controller will give an audible signal!

Set the controller power switch to the OFF position.

Plug the controller power cord into a 115V, 50/60 Hz or 230 V, 50/60 Hz grounded AC power supply, as required.

SYSTEM ALIGNMENT & SPINDLE INSERTION

Looking down on the viscometer, level the instrument by adjusting the leveling screws on the base until the viscometers bubble level is centered.

Position the Thermo Container and base between the viscometer stand legs and adjust the three (3) knurled screws until the bubble level is centered on the thermosel.

Lower the viscometer until the tips of the alignment bracket (see Figure I-1) just touch the horizontal surface of the locating ring, making contact directly behind the vertical

curve. Raise the viscometer, positioning the tips of the alignment bracket about 1/16 inch above the horizontal surface behind the locating ring. An etched line on the back of the locating ring is the 1/16 inch reference point.

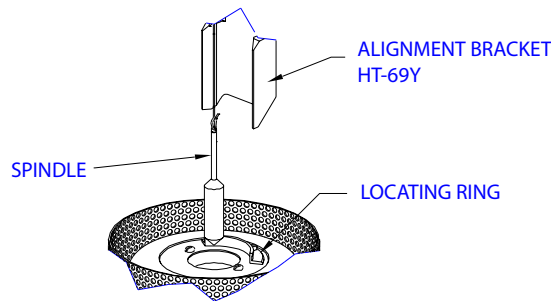


Figure I-1

NOTE: Do not forcibly displace the alignment bracket!

Add the appropriate sample volume to the chamber (see Appendix A, Range Tables, for volume requirement; see viscometer operating manual for calibration procedures). Carefully pour, pipette, or syringe the sample into the chamber; avoid introducing air bubbles; tilt the chamber slightly by hand to facilitate the pouring process. Place the chamber into the Thermo Container.

- If using a spindle that requires an extension link and coupling nut, attach one end of the link to the spindle and the other to the coupling nut. Carefully lower the spindle by the coupling nut and link into the sample chamber. While pushing up gently on the pivot screw that protrudes from the bottom of the viscometer, thread the coupling nut onto the viscometer. Note that the spindle has a left-hand thread.
- If using a Magnetic spindle, attach the spindle coupling to the viscometer by carefully inserting the spindle coupling into the instrument coupling. There should be no gap between the spindle coupling and the instrument coupling, when properly installed.
- If using a solid shaft (SC4-**BS) spindle, carefully lower the spindle into the sample chamber. Install the spindle coupling to the instrument as noted above.

Place the insulating cap over the sample chamber inlet.

Brookfield Thermosel with DV2T Viscometer

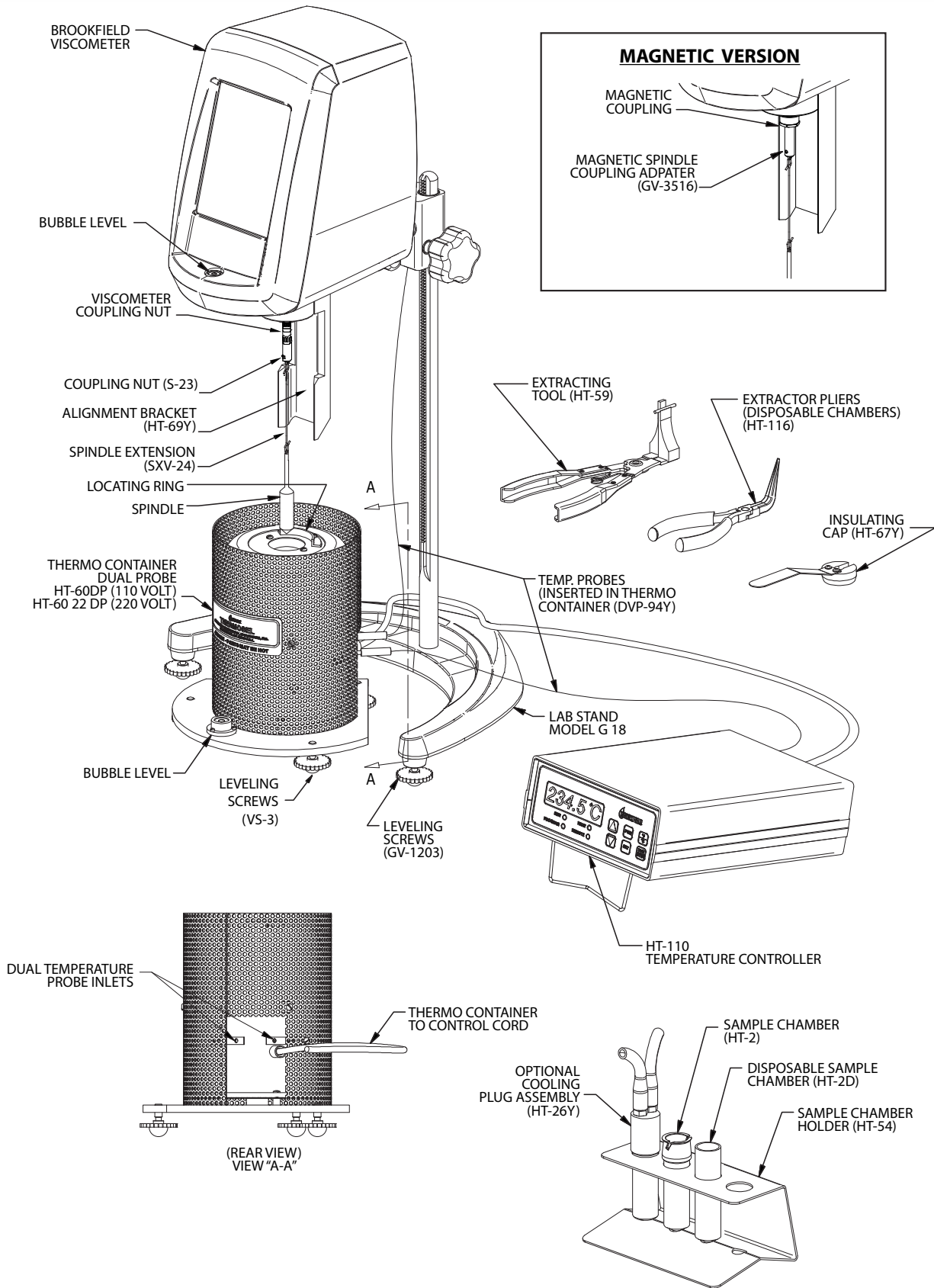


Figure I-2

I.5 GENERAL

The operator should become familiar with the alignment procedure so that the system can be safely aligned and operated at elevated temperatures.

The system is designed to operate in the temperature range from 40°C (104°F) to a maximum of 300°C (572°F). Minimum test temperature when using the ThermoSel should be 20°C above ambient temperature. Control of test sample temperature is possible in this range with accuracy as follows:

Temperature Range	Temperature Accuracy
40°C to +150°C	±1.0°C
+150°C to +300°C	±2.0°C

NOTE: *The lower limit of control is approximately 40°C (104°F). The upper limit of precise control is 300°C (572°F). This is determined not by the controller-heat element capabilities, but by design of the spindle, sample chamber, thermo-container, and the proximity of the viscometer to the thermo-container.*



Use of the system above 300°C could result in damage to the Thermo Container and to the viscometer.

The controller will bring the Thermo Container to elevated set point temperature in approximately 30 minutes. However, lowering the temperature from the elevated set point can be quite time consuming. Approximately four hours is required for the Thermo Container to cool from 260°C (500°F) to 37°C (100°F). Using the optional cooling plug (HT-26Y), this temperature reduction can be accomplished in about 20 minutes (see Figure I-3). Familiarization with the temperature controller instructions should be made at this time.

Refer to Appendix A for viscosity range data, spindle factors (for use with Dial Reading Viscometers), and test volume sample size. This information is to be used when obtaining viscosity measurements.

I.6 VISCOSITY MEASUREMENTS

SAMPLES THAT ARE LIQUID AT AMBIENT TEMPERATURE

With the instrumentation assembled and familiarization completed, the following steps are recommended for taking viscosity measurements of samples which are liquid at ambient temperature:

1. Remove the insulating cap; for a standard spindle, unthread the coupling nut and remove the spindle; for a Magnetic spindle, lift the spindle up and push the spindle to the side to disengage it.
2. The solid shaft magnetic spindle can be disengaged by raising the spindle out of the ThermoSel and then pushing the spindle to the side to disengage it.
3. Raise the viscometer to the highest level on the stand.
4. Remove the sample chamber (HT-2 or HT-2D) using the extracting tool (HT-59 or

HT-116 respectively) and place it in the sample chamber holder (HT-54).

5. Using a syringe, graduated cylinder, or other suitable measuring device, pour into the sample chamber the volume of liquid sample specified on the range table in Appendix A.

NOTE: *Avoid introducing air bubbles into the sample.*

6. Using an extracting tool, put the filled sample chamber back into the Thermo Container by rotating the chamber until it drops and locks in place.
7. Lower the viscometer and align the thermo-container.
8. Lower the spindle into the chamber and connect the link and coupling nut (S-23) to viscometer by carefully lifting the viscometer shaft up slightly while screwing on the S-23. (Note: left-hand thread). Or for a Magnetic version instrument, connect the spindle coupling (GV-3516) to the instrument by carefully inserting the spindle coupling into the instrument coupling. There should be no gap between the spindle coupling and the instrument coupling, when properly installed.
9. Check that the liquid level is approximately 1/8 inch above the conical surface on the spindle shaft. Do not overfill.
10. Replace the insulating cap.
11. Turn on the temperature controller and adjust the set point to the desired temperature.

NOTE: *The set point temperature will be denoted by a steady display on the controller, along with the steady flashing of the heater “ON” light.*

One option for sample materials that are not thixotropic or pseudoplastic is to turn ON the viscometer motor and leave it running during the equilibrium period. The rotating spindle provides agitation which helps reduce temperature equilibrium time and temperature gradients within the test sample.

12. After the Thermo Container, spindle, chamber, and test sample have reached temperature equilibrium, viscosity readings may be obtained.

SAMPLES THAT ARE NOT LIQUID AT AMBIENT TEMPERATURE

There are several alternative approaches toward testing samples that are solid or semi-solid at ambient (room) temperature. The main concerns are: First - safety; Second - keeping test material IN the sample chamber and OUT OF the Thermo Container “well.” The Thermo Container can be preheated without the sample chamber and test material if desired.

The following steps are recommended for taking viscosity measurements of test samples that are liquid only at elevated temperatures.

1. The instrumentation should be assembled without the insertion of the sample chamber, spindle, and insulator cap.
2. Set the controller to the desired temperature.

3. Weigh into the sample chamber the amount of test sample which in the liquid state will be equivalent to the test volume required. It is helpful to know the density of the sample material.
4. Place the sample chamber into the Thermo Container and proceed, following the steps outlined for liquid samples at ambient temperature, starting on Page 6.

THERMO CONTAINER ASSISTED COOLING

With the sample chamber removed, the Thermo Container can be cooled by inserting the optional cooling plug (HT-26Y) into the sample chamber well and circulating a cooling medium (tap water) through it.



Water must be circulating through the cooling plug at all times prior to and during the cooling period.

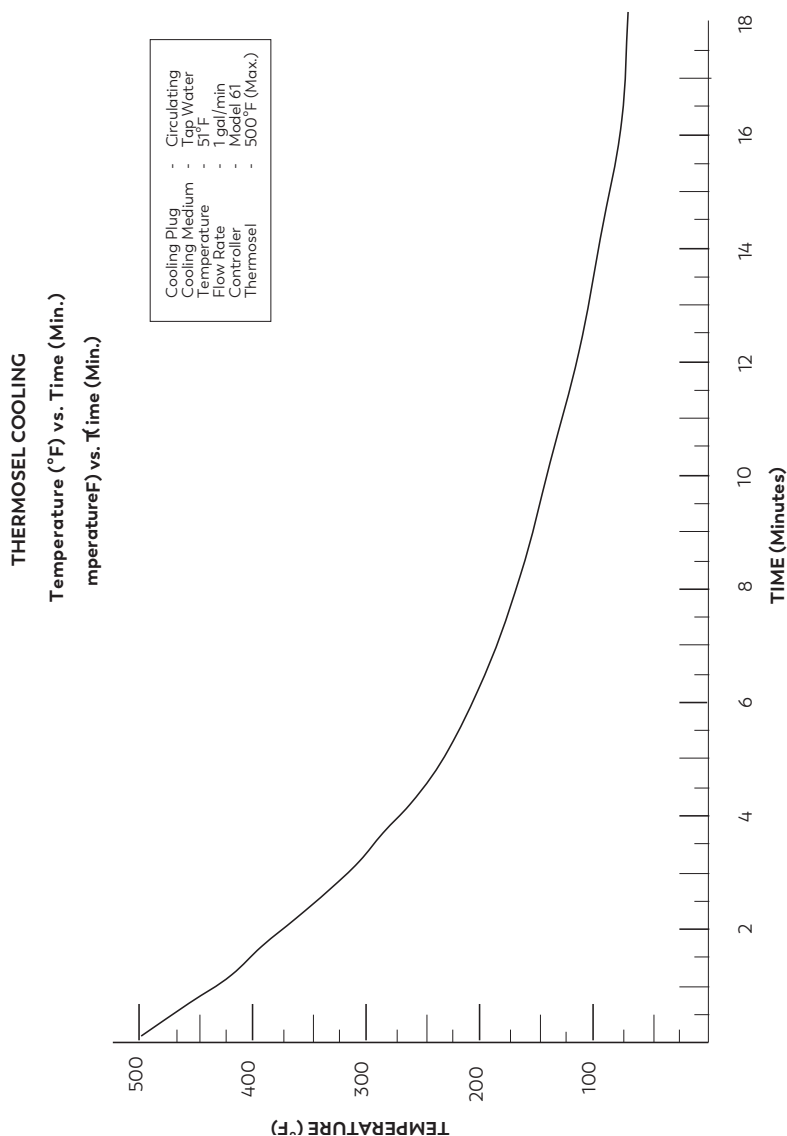


Figure I-3

I.7 CLEAN-UP

CLEANING

Using the extracting tool, lift the chamber out of the Thermo Container and pour out the sample material. The spindle and chamber are made of stainless steel which can be cleaned with most commercial solvents. They should not come in contact with Sulfuric Acid; Hydrofluoric Acid; Ferric Chloride or Hydrochloric Acid solutions. Use mechanical action as required to scrape out the chamber.

Disposable sample chambers and spindles do not require any type of cleaning and may be discarded after use.

Regulated materials that are hazardous require proper disposal in accordance with local regulations.

APPENDIX A - VISCOSITY RANGES / SPINDLE FACTORS

THERMOSEL VISCOSITY RANGE DATA

DIAL VISCOMETERS

Spindle	Viscosity (cP)		Shear Rate sec ⁻¹	Sample Volume (mL)
	LVT			
SC4-18	5	- 10,000	1.32N	8.0
SC4-31	50	- 100,000	0.34N	10
SC4-34	100	- 200,000	0.28N	9.5
HT-DIN-81	5.7	- 11,400	1.29N	7.0

Spindle	Viscosity (cP)		Shear Rate sec ⁻¹	Sample Volume (mL)
	RVT			
SC4-21	50	- 100,000	0.93N	8.0
SC4-27	250	- 500,000	0.34N	10.5
SC4-28	500	- 1,000,000	0.28N	11.5
SC4-29	1,000	- 2,000,000	0.25N	13.0
HT-DIN-81	36.5	- 73,000	1.29N	7.0

Spindle	Viscosity (cP)		Shear Rate sec ⁻¹	Sample Volume (mL)
	HAT	HBT		
SC4-21	100 - 200,000	400 - 800,000	0.93N	8.0
SC4-27	500 - 1,000,000	2,000 - 4,000,000	0.34N	10.5
SC4-28	1,000 - 2,000,000	4,000 - 8,000,000	0.28N	11.5
SC4-29	2,000 - 4,000,000	8,000 - 16,000,000	0.25N	13.0
HT-DIN-81	73 - 146,000	292 - 584,000	1.29N	7.0

N = RPM

DVE, DV-E, DV1 AND DV-I SERIES VISCOMETERS

Spindle	Viscosity (cP)	Shear Rate sec ⁻¹	Sample Volume (mL)
	LV Torque		
SC4-18	3 - 10,000	1.32N	8.0
SC4-31	30 - 100,000	0.34N	10
SC4-34	60 - 200,000	0.28N	9.5
HT-DIN-81	3.42 - 11,400	1.29N	7.0

Spindle	Viscosity (cP)			Shear Rate sec ⁻¹	Sample Volume (mL)
	RV Torque	HA Torque	HB Torque		
SC4-21	50 - 170,000	100 - 334,000	400 - 1,336,000	0.93N	8.0
SC4-27	250 - 830,000	500 - 1,660,000	2,000 - 6,640,000	0.34N	10.5
SC4-28	500 - 1,660,000	1,000 - 3,320,000	4,000 - 13,280,000	0.28N	11.5
SC4-29	1,000 - 3,330,000	2,000 - 6,660,000	8,000 - 26,640,000	0.25N	13.0
HT-DIN-81	36.5 - 121,700	73 - 243,300	292 - 973,300	1.29N	7.0

N = RPM

DV2T AND DV-II SERIES VISCOMETERS

Spindle	Viscosity (cP)	Shear Rate sec ⁻¹	Sample Volume (mL)
	LV Torque		
SC4-18	3 - 30,000	1.32N	8.0
SC4-31	15 - 300,000	0.34N	10
SC4-34	30 - 600,000	0.28N	9.5
HT-DIN-81	3.42 - 34,200	1.29N	7.0

Spindle	Viscosity (cP)			Shear Rate sec ⁻¹	Sample Volume (mL)
	RV Torque	HA Torque	HB Torque		
SC4-21	25 - 500,000	50 - 1,000,000	200 - 4,000,000	0.93N	8.0
SC4-27	125 - 2,500,000	250 - 5,000,000	1,000 - 20,000,000	0.34N	10.5
SC4-28	250 - 5,000,000	500 - 10,000,000	2,000 - 40,000,000	0.28N	11.5
SC4-29	500 - 10,000,000	1,000 - 20,000,000	4,000 - 80,000,000	0.25N	13.0
HT-DIN-81	18.25 - 365,000	36.5 - 730,000	146 - 2,920,000	1.29N	7.0

N = RPM

DV3T & DV-III SERIES RHEOMETERS

Spindle	Viscosity (cP)	Shear Rate sec ⁻¹	Sample Volume (mL)
	LV Torque		
SC4-18	3 - 30,000	1.32N	8.0
SC4-31	12 - 300,000	0.34N	10
SC4-34	24 - 600,000	0.28N	9.5
HT-DIN-81	3.42 - 34,200	1.29N	7.0

Spindle	Viscosity (cP)			Shear Rate sec ⁻¹	Sample Volume (mL)
	RV Torque	HA Torque	HB Torque		
SC4-21	20 - 500,000	40 - 1,000,000	160 - 4,000,000	0.93N	8.0
SC4-27	100 - 2,500,000	200 - 5,000,000	800 - 20,000,000	0.34N	10.5
SC4-28	200 - 5,000,000	400 - 10,000,000	1,600 - 40,000,000	0.28N	11.5
SC4-29	400 - 10,000,000	800 - 20,000,000	3,200 - 80,000,000	0.25N	13.0
HT-DIN-81	14.6 - 365,000	29.2 - 730,000	117 - 2,920,000	1.29N	7.0

N = RPM

THERMOSEL SPINDLE FACTORS

SPEED (RPM)	LV VISCOMETER SPINDLE NUMBER		
	18	31	34
60	0.5	5	10
30	1	10	20
12	2.5	25	50
6	5	50	100
3	10	100	200
1.5	20	200	400
0.6	50	500	1K
0.3	100	1K	2K

SPEED (RPM)	RV VISCOMETER SPINDLE NUMBER			
	21	27	28	29
100	5	25	50	100
50	10	50	100	200
20	25	125	250	500
10	50	250	500	1K
5	100	500	1K	2K
4	125	625	1.25K	2.5K
2.5	200	1K	2K	4K
2	250	1.25K	2.5K	5K
1	500	2.5K	5K	10K
0.5	1K	5K	10K	20K

SPEED (RPM)	HA VISCOMETER SPINDLE NUMBER			
	21	27	28	29
100	10	50	100	200
50	20	100	200	400
20	50	250	500	1K
10	100	500	1K	2K
5	200	1K	2K	4K
2.5	400	2K	4K	8K
2	500	2.5K	5K	10K
1	1K	5K	10K	20K
0.5	2K	10K	20K	40K

SPEED (RPM)	HB VISCOMETER SPINDLE NUMBER			
	21	27	28	29
100	40	200	400	800
50	80	400	800	1.6K
20	200	1K	2K	4K
10	400	2K	4K	8K
5	800	4K	8K	16K
2.5	1.6K	8K	16K	32K
2	2K	10K	20K	40K
1	4K	20K	40K	80K
0.5	8K	40K	80K	160K

K=1000 To calculate viscosity in centipoise (cP), multiply the dial reading by the factor corresponding to the viscometer spindle and speed combination utilized.

APPENDIX B - HIGH TEMPERATURE CALIBRATION FLUIDS

These fluids are recommended for verifying the calibration of the Brookfield Thermosel System.

Fluid	Temperature		Nominal Viscosity Centipoise (mPa•s)
	°C	°F	
HT - 30,000	25.0	77	30,000
	93.3	200	9,000
	149.0	300	4,500
HT-60,000	25.0	77	60,000
	93.3	200	18,000
	149.0	300	9,000
HT-100,000	25.0	77	100,000
	93.3	200	30,000
	149.0	300	15,000

The viscometer calibration is verified first. A regular* “disc” spindle should be used to test the HT fluid in its jar at 25.0°C or 77°F. Please refer to the appropriate viscometer operation manual for instructions.

The Thermosel System and its appropriate spindle are then used to measure the HT fluid’s viscosity at 200 and 300°F.

Contact AMETEK Brookfield or an authorized dealer in your area for purchase of these fluids.

* Standard spindles supplied with viscometer (not RV/HB-7 or LV-4 spindles).

APPENDIX C - DISPOSABLE SAMPLE CHAMBER AND SPINDLES

Disposable Sample Chambers are intended for use with difficult to clean materials, or in busy laboratories where the low cost disposable chamber may improve productivity. Also available are disposable SC4-27D spindles for use with the Brookfield Thermosel.

The HT-2DB disposable chambers are inserted into or removed from HT-60DP or HT-60ADP Thermo Container using the HT-116 pliers as shown in Figure C-1.

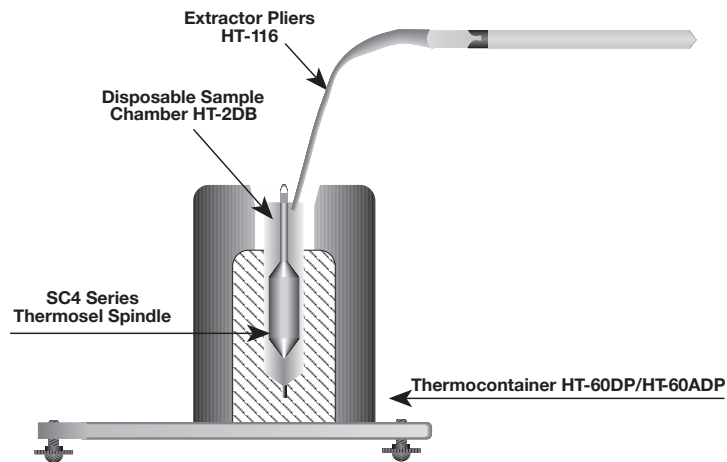


Figure C-1

Replacement disposable chambers are available in packages of 100 chambers (Brookfield Part No. HT-2DB-100). Contact AMETEK Brookfield or your local authorized dealer for pricing and delivery information.

A package of five disposable aluminum sample chambers (Part No. HT-2DB) and one pair of extraction pliers (Part No. HT-116) are supplied with each new Thermosel System.

DISPOSABLE SPINDLE INSTRUCTIONS (SEE FIGURES C-2 AND C-3)

1. Loosen closer.
2. Slide disposable spindle into chuck to insertion mark.
3. Hold chuck and hand-tighten closer.
4. For standard assembly, attach the assembly (spindle and chuck) to the viscometer being careful to hold the coupling nut firmly and lift up slightly while screwing on the assembly (Note: left-handed thread). See Figure C-2.
5. For Magnetic assembly, attach the assembly (spindle and chuck) to the viscometer by carefully inserting the spindle coupling into the instrument coupling. There should be no gap between the spindle coupling and the instrument coupling, when properly installed. See Figure C-3.

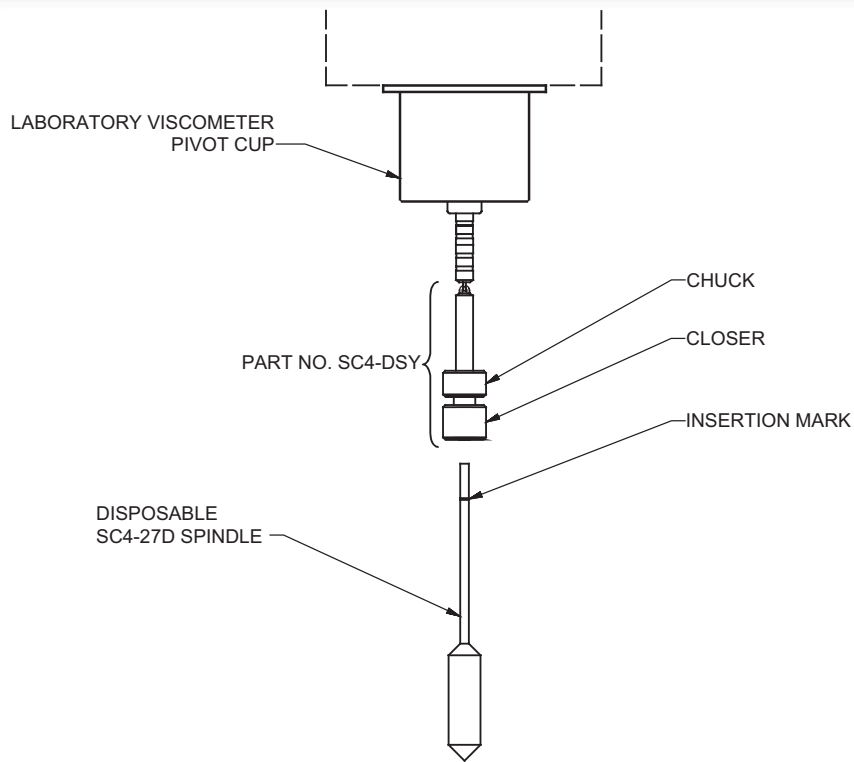


Figure C-2: Standard

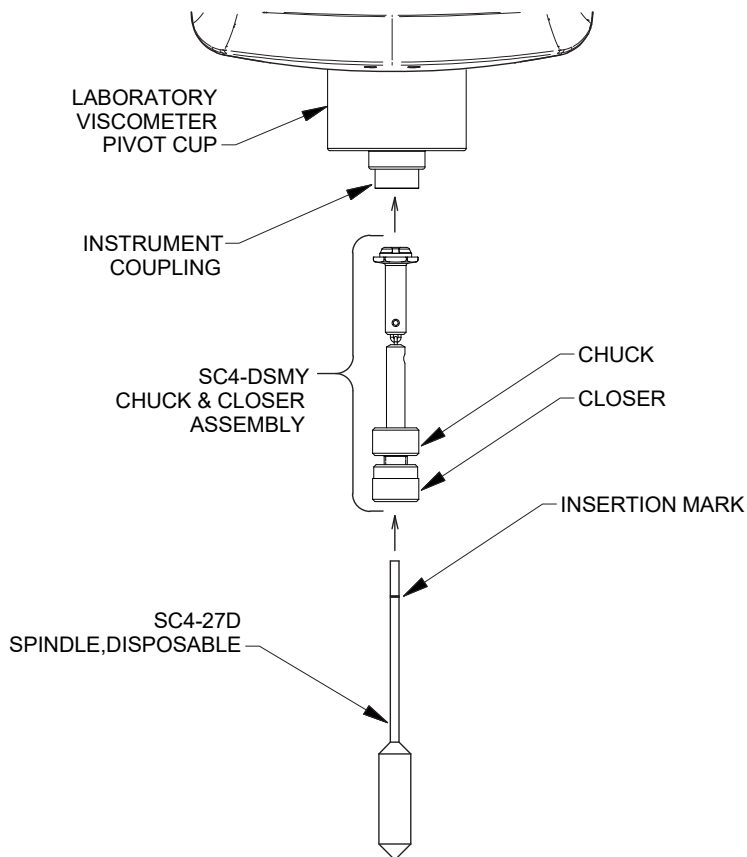


Figure C-3: Magnetic

APPENDIX D - SOLID SHAFT SPINDLE OPTION

Solid shaft spindles are available for use with the Brookfield Thermosel Systems. These spindles facilitate the measurement of “stiff”, slow-flowing materials such as:

Candy base
Highly-filled epoxies

Pastes
Asphalts

The solid shaft on these spindles minimizes the “walking” (eccentric rotation) experienced when measuring “stiff” materials with the standard link hanging Thermosel spindles. These spindles are directly interchangeable in viscosity range and sample volume with the standard Thermosel spindles. No modifications to the Thermosel or the viscometer are necessary. Solid shaft spindles are available, as listed:

Standard Solid Shaft Spindles	Magnetic Solid Shaft Spindles
SC4-27 BS	SC4-27 MBS
SC4-28 BS	SC4-28 MBS
SC4-29 BS	SC4-29 MBS
SC4-31 BS	SC4-31 MBS
SC4-34 BS	SC4-34 MBS

Contact AMETEK Brookfield or your local authorized dealer for purchase of these spindles.

APPENDIX E - ONLINE HELP AND ADDITIONAL RESOURCES

[WWW.BROOKFIELDENGINEERING.COM](http://www.brookfieldengineering.com)

The Brookfield website is a good resource for additional and self-help whenever you need it. Our website offers a selection of “how-to” videos, application notes, conversion tables, instructional manuals, material safety data sheets, calibration templates and other technical resources.

[HTTP://WWW.YOUTUBE.COM/USER/BROOKFIELDENG](http://www.youtube.com/user/BROOKFIELDENG)

Brookfield has its own YouTube channel. Videos posted to our website can be found here as well as other “home-made” videos made by our own technical sales group.

MORE SOLUTIONS TO STICKY PROBLEMS

Learn more about viscosity and rheology with our most popular publication. This informative booklet will provide you with measurement techniques, advice and much more. It’s a must-have for any Brookfield Viscometer or Rheometer operator. More Solutions is available as a downloadable PDF on the Brookfield website by following this path: <https://www.brookfield-engineering.com/downloads/technical-documents>.

TRAINING/COURSES

Whether it is instrument-specific courses, training to help you better prepare for auditing concerns, or just a better understanding of your methods, who better to learn from than the worldwide leaders of viscosity measuring equipment? Visit our [Services](#) section on our website to learn more about training.

APPENDIX F - WARRANTY REPAIR AND SERVICE

Brookfield Viscometers/Rheometers are guaranteed for one year from date of purchase against defects in materials and workmanship. They are certified against primary viscosity standards traceable to the National Institute of Standards and Technology (N.I.S.T.). The Rheometer must be returned to AMETEK Brookfield or to the authorized dealer from whom it was purchased for a warranty evaluation. Transportation is at the purchaser's expense. The Rheometer should be shipped in its carrying case together with all spindles originally provided with the instrument. If returning to Brookfield, please contact us for a return authorization number prior to shipping.

All Brookfield DVNext Rheometers are supplied from the factory with a Calibration Seal (located on the back of the viscometer). The warranty stated above will be voided if the Calibration Seal has been damaged. Only Brookfield or our authorized servicing dealers may break the Calibration Seal for purposes of instrument repair or recalibration.

For a copy of the Repair Return Form, go to the Brookfield website,
<https://www.brookfieldengineering.com/>

For repair or service in the United States return to:

AMETEK Brookfield
11 Commerce Boulevard
Middleboro, MA 02346 U.S.A.
Telephone: (508) 946-6200 FAX: (508) 946-6262
www.brookfieldengineering.com

For repair or service outside the United States, consult AMETEK Brookfield or the local authorized dealer from whom you purchased the instrument.

For repair or service in the United Kingdom return to:

AMETEK (GB) Limited
Brookfield Technical Centre
Stadium Way
Harlow, Essex CM19 5GX, England
Telephone: (44) 1279/451774 FAX: (44) 1279/451775
www.brookfieldengineering.uk

For repair or service in Germany return to:

AMETEK GmbH
Oberweyerer Straße 21
65589 Hadamar-Steinbach,
Telephone: (49) 7172/927100 FAX: (49) 7172/927105
www.brookfieldengineering.de

For repair or service in China return to:

AMETEK Commercial Enterprise (Shanghai) Co., Ltd Guangzhou Branch
Room 810 Dongbao Plaza, No. 767 East Dongfend Road
Guangzhou, 510600 China
Telephone: (86) 20/3760-0548 FAX: (86) 20/3760-0548
www.brookfieldengineering.cn

On-site service at your facility is also available from Brookfield. Please contact our Service Department in the United States, United Kingdom, Germany or China for details.